

Instructions for Faucet Mounted Eye Wash Devices Models: 0910, 0920, and 0930

Installation

1. Remove the aerator or other fitting from faucet outlet.
2. Mount the eye wash unit to faucet spout. The eye wash inlet has a 3/4" x 27 female thread.

Note: Due to the wide variety of faucets, it may be necessary to use an adapter to mount the eye wash unit on the faucet spout. Three adapters are furnished with the unit: the 400-10 (larger) adapter has a 15/16" X 27 male thread to fit Delta and American Standard faucets, the 400-9 (smaller) adapter has a 13/16" x 27 thread to fit Kohler and other faucets, and the 400-11 adapter has a 3/8" IPS male thread to fit laboratory type faucets. Additional adapters may be purchased at most hardware stores.

3. On models 0910 and 0920, adjust the angle of the outlet heads to obtain the desired angle for the eye wash water flow by loosening the outlet arm nut, positioning the heads as desired by re-tightening the nut. If an insufficient flow of water is delivered by the unit, the flow may be increased by removing one or more screens from the eye wash outlet heads aerators.
4. Attach the enclosed sign to wall or backsplash behind sink.

Use

1. Turn on the cold water side of the faucet. **Important: use cold water only when the eye wash unit is in operation. Use of hot water may cause scalding of the eye tissue.**
2. Pull the knob on the eye wash unit forward to activate the eye wash. Water pressure will hold the eye wash in operation, leaving the user's hands free.
3. To return to normal faucet operation, push knob back to the original position. When the faucet is turned off, the knob will return to the original position automatically.

Maintenance

1. Periodically clean the eye wash aerators and valve to assure proper water flow.
2. Keep plastic float-off covers on outlet heads when the unit is not in use.
3. The eye wash unit, like all emergency eye wash and shower equipment, should be tested weekly.

Note: Faucet mounted eye wash units require two motions to operate (turn faucet on, pull eye wash knob). Because they cannot be activated in one second or less, these units do not comply with ANSI Z358.1-1990. Therefore, faucet mounted eye wash units are intended as supplements to, but not replacements for, standard plumbed or self contained eye wash.

fend·all
World leaders in emergency eyecare.

5 East College Drive, Arlington Heights, IL 60004
847-577-7400 • 800-543-4842